

*Spare parts list*  
*BIOBASIC RF 310*




Description :

Spare parts list for  
 BIOBASIC RF 310

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
03.09.15	JP	30.09.22	JP	008		

Exploded view

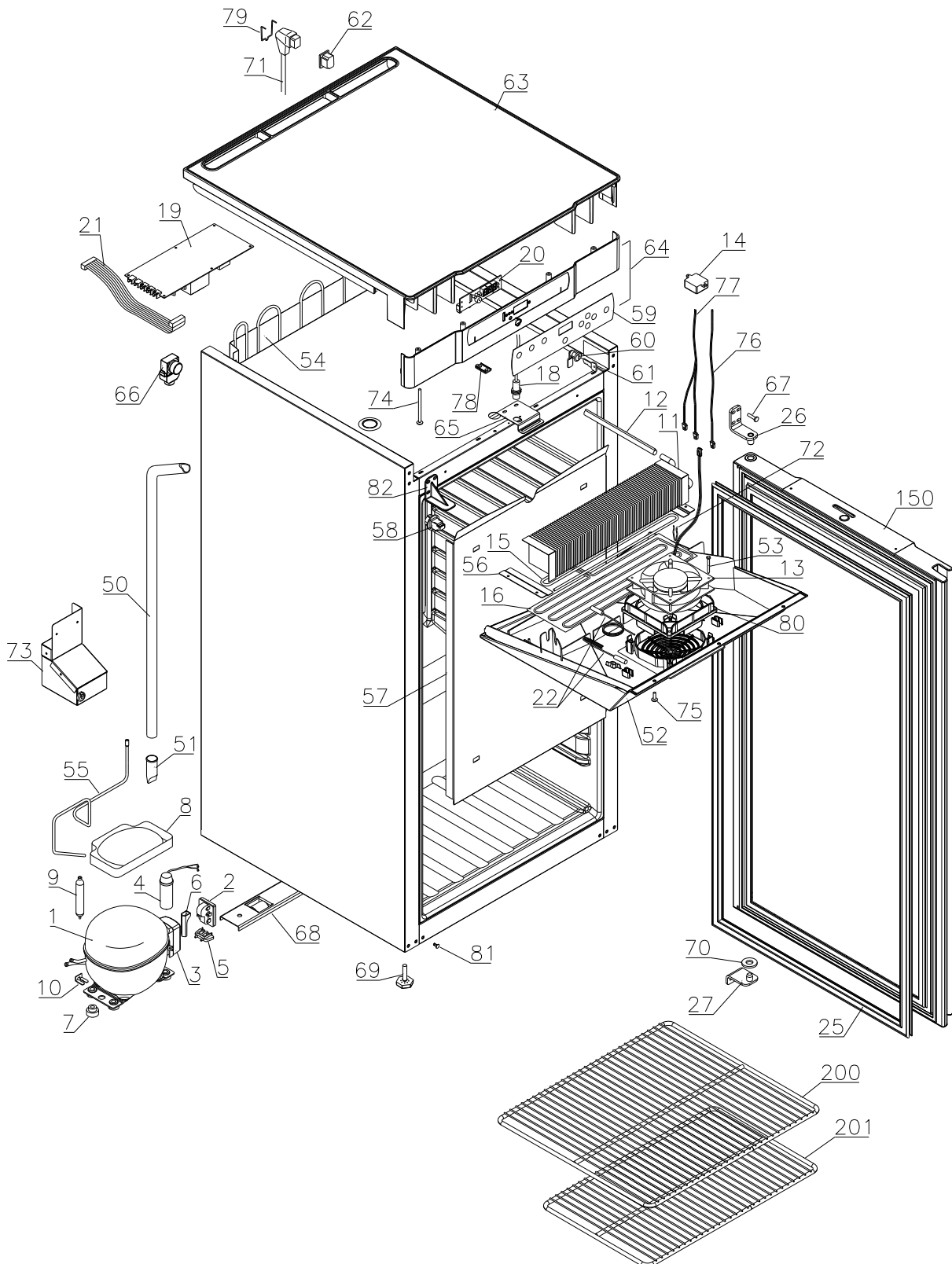

**Gram Commercial**  
**Technical information**

765041763

Description :

**Spare parts list for  
BIOBASIC RF 310**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
03.09.15	JP	30.09.22	JP	008		

Item no.	Description	Gram part no.	Week/year:	
			Issued	Deleted

**Essential spare parts**

RF 310 G

1	Compressor NLX13KK.1	500550096	3715	
2	Starting device	453050178	3715	
3	Cover	452440293	3715	
4	Run capacitor 4 µF	451723513	3715	
5	Cord relief	452440297	3715	
6	Protection screen PTC	452440290	3715	

RF 310 H

1	Compressor NLY9FK	500550095	3715	
2	Starting device	453050178	3715	
3	Cover	452440293	3715	
4	Run capacitor 4µF	451723513	3715	
5	Cord relief	452440297	3715	
6	Protection screen PTC	452440290	3715	

Common parts:

7	Rubber bush	760080519	3715	
8	Re-evaporation tray	410820011	3715	
9	Filter drier	761221322	3715	
10	Compressor clip	760082367	3715	
11	Evaporator	765080207	3715	
12	Heat exchanger	761221406	3715	
13	Evaporator fan 12V DC	765071196	3715	
14	Trafo for fan (kit)	761234127	3715	1521
14	Trafo for fan	765072987	1521	
15	Defrost heater	765071197	3715	
16	Drip tray heater	765071199	3715	
18	Door switch	450400111	3715	
19	Electr.contr., power part F70	761256165	3715	
20	Electronic controller, display part	612900056	3715	
21	Band cable	765072661	3715	
22	Temperature sensor	453531501	3715	
25	Door gasket	761222728	3715	
26	Top hinge	761222906	3715	
27	Bottom hinge	761222904	3715	

**Gram Commercial**  
**Technical information**

765041763

Description :

**Spare parts list for  
BIOBASIC RF 310**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
03.09.15	JP	30.09.22	JP	008		

Item no.	Description	Gram part no.	Week/year Issued	Deleted
----------	-------------	---------------	---------------------	---------

**Other spare parts**

50	Drain pipe	761221758	3715	
51	Water trap	760082891	3715	
52	Drip tray, white	761223101	3715	
53	Screw for fan	420790001	3715	
54	Condenser	761222786	3715	
55	Re-evaporation pipework	760000323	3715	
56	Evaporator suspension bracket	761223070	3715	
57	Air distribution plate	761222994	3715	
58	Suspension plug, white	765081213	3715	
59	Label for controller	765041713	3715	
60	Door lock	765080093	3715	
61	Key	761220166	3715	
62	Appliance inlet	451022002	3715	
63	Tabletop upper part, white	761222630	3715	
64	Tabletop front, white	761222414	3715	
65	Lock bracket	761222761	3715	
66	Air relief valve	765081206	3715	
67	Screw for hinge	420090100	3715	
68	Wheel bracket complete	761223051	3715	
69	Adjustment screw	761223055	3715	
70	Washer	760250524	3715	
71	Mains cable, EU	765072630	3715	
71	Mains cable, UK	765072631	3715	
71	Mains cable, DK	765072632	3715	
72	Bracket for evaporator heater	761222779	3715	
73	Condenser fan complete	761221780	3715	

*Gram Commercial*  
*Technical information*

765041763

Description :

**Spare parts list for  
BIOBASIC RF 310**

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
03.09.15	JP	30.09.22	JP	008		

Item no.	Description	Gram part no.	Week/year Issued	Deleted
74	Screw M4x70	420178125	3715	
75	Screw M4x16	422701019	3715	
76	Fan wire	765072656	3715	
77	Heater wire	765073327	3715	
78	Cover plug, white	765081247	3715	
79	Brace for mains cable	761230215	3715	
80	Rubber gasket	761223110	3715	
81	Cover plug, white	765080504	3715	
82	Cover bracket, white	761222758	3715	
150	Door complete, white	761232759	3715	
200	Wire shelf large, white	818721001	3715	
201	Wire shelf small, white	818721019	3715	